LED-UV-Insektenleuchte
Entomology LED UV Lamp

entoLED[®] mini

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Vielen Dank

General Information

This high quality product has been designed and developed to require a minimum of maintenance. Please follow the guidelines below and read this manual in its entirety to ensure safe operation of the unit.

Symbols

The following symbols are used in this manual.

CAUTION! This symbol indicates a potential risk and alerts you to proceed with caution.

CAUTION! This symbol indicates a potential danger of optical radiation.

Warning

- Please always disconnect mains plug before cleaning the device.

Safety Information

- To avoid eye damage, don't look directly into the light source when turned on.
- Versions with UV light release intensive ultraviolet radiation during operation.

A major part of the light is invisible. The beams can damage your health. Protect your eyes (UV protection glasses) and skin during operation.

Important: The evaluation of eye safety occurs according to the standard CIE S009/E:2002 („photo biological safety of lamps and lamp systems“). Within the risk grouping system of this CIE standard, the LEDs specified in this data sheet fall into the „low risk“ group (relating to devices in the visible spectrum with an exposure time of 100s). Under real circumstances (for exposure time, eye pupils, observation distance), it can be assumed that no endangerment to the eye exists from these devices. As a matter of principle, however, it should be mentioned that intense light sources have a high secondary exposure potential due to their blinding effect. As is also true when viewing other bright light sources (e.g. headlights), temporary reduction in visual acuity and afterimages can occur, leading to irritation, annoyance, visual impairment, and even accidents, depending on the situation.

Function and Operation

Package Contents

Entomology lamp entoLED mini	1
Multi adapter power supply (12 V, EU UK US)	1
Connection cable with battery clamps	1
Protective travel pouch	1
User Manual	1

Technical Data

LED Type	5 High Power LEDs
Illuminants	2 × UV (365 nm), 1 × UV (375 nm), 1 × UV (385 nm), 1 × royal blue (450 nm)
Diameter	64 mm
Height	30 mm
IP Code	IP65
Reverse polarity protection	yes
Power consumption	approx. 3,7 W
Operating voltage with power supply	100 - 240 V AC, 50/60 Hz
Operating voltage entoLED	12 V DC
Color	olive green
Material	anodized aluminium, borosilicate glass
Weight	approx. 240 g

Getting Started


The LED lamp is a plug & play device very easy to use.

Please check before use whether...

...the mains plug is connected.

...the environment light is dark enough for the optionally day light sensor, cover these if necessary.

Description of Buttons and Switches

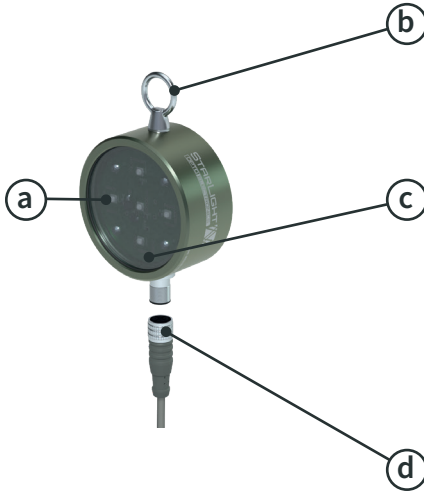- a) **Illuminants**
- b) **Mounting lug**
- c) **Borosilicate glass**
- d) **Connection from the power supply or battery**
Red +
Black -

Technical Assistance

Troubleshooting

The information in the table below can help to resolve the possible operating problems quickly.

Problem	Causes	Solution
No function at the device	no power connection	Please check whether the power cables are connected.
	electronic defective	Please contact your distributor.
	LED defective	Please contact your distributor.
LED gets dark	thermal overload of the system	Please ensure an ambient temperature from 5° to 35°C.
	LED life span exceeded	Please contact your distributor.

Storage, Maintenance, Repair

The device is maintenance-free. Some points, however, you should take into consideration.

1. Please unplug the device before cleaning.
2. For the cleaning of the housing please use a smooth, semi-wet cloth, if necessary with a smooth household cleaning fluid.
3. No moisture should entering into the housing interior.
4. To avoid damages during storage or transport please retain the packaging.

Repairs are only allowed by trained and experienced personnel. In case of improper handling the warranty expires.

Technical Service

Web Resources

Visit our website at entoLED.eu for:

- Contact details of technical customer service.
- Product overview as well as infos about accessories and spare parts.
- Additional product information and special offers.

Contact

For more information or technical assistance please contact:

bioform entomology

Dr. J. Schmidl e.K.

Am Kressenstein 48 90427 Nürnberg (Germany)

www.entoLED.eu

info@bioform.de

Warranty

StarLight Opto-Electronics GmbH & Co. KG warrants that this product will be free from defects in material and workmanship for a period of five (5) years from date of purchase. If a defect is present, StarLight will, at its option, repair, replace, or refund the purchase price of this product at no charge to you. Precondition is the device is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, or misapplication, or from ordinary wear and tear.

For your protection, items being returned must be insured against possible damage or loss. This warranty shall be limited to the replacement of defective products.

It is expressly agreed that this warranty will be in lieu of all warranties of fitness and in lieu of the warranty of merchantability.

Equipment Disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste. Instead, it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you

Notizen / Notes